



Certified Balance & Scale

www.certify911.com

office@certify911.com

ph: 800-777-1198

ph: 815-895-5158

Quick start guide: YSI pH 1110 meter

Do the following to set up:

1. Plug in probe to back of meter. (Temp sensor will plug into top hole "TP")
2. Plug adapter into back of meter and plug in to outlet

To use meter:

Follow the calibration instructions in manual or calibration guide book (multi-language book). Here is a quick walk through:

Turn on meter

To calibrate

1. Press [Cal] "ct 1" displays
2. Rinse probe and place in first buffer (usually pH 7).
3. Wait 30 seconds. Press [Enter]

After meter calibrates 7.0 it will display "ct2".

4. Rinse probe with DI water and then place in second buffer.
If you are doing a 2 pt., use the pH that is in your range (4 or 10).
5. Wait 30 seconds. Press [Enter]

After meter calibrates it will display "ct3".

6. If doing a 2 pt. Calibration, press [M] and then [Enter] to display the slope.

To read a sample, press [M] to return to pH measurement mode.

Always store the probe in the solution of 3mol KCL solution. If not available, use pH 4.0

Contact us for supplies, equipment, and yearly certification options:

www.certify911.com