



Certified Balance & Scale

www.certify911.com

office@certify911.com

ph: 800-777-1198

ph: 815-895-5158

Quick start guide: YSI pH 1110 meter

Do the following to set up:

1. Plug in probe to back of meter. (Temp sensor will plug into top hole "TP")
2. Plug adapter into back of meter and plug in to outlet

To use meter:

Follow the calibration instructions in manual or calibration guide book (multi-language book). Here is a quick walk through:

1. Turn on meter
2. Press [Cal] until "AutoCal TEC" displays
3. Rinse probe with buffer 7.0 and place in buffer.
4. Wait 30 seconds. Press [Enter]
5. When meter calibrates 7.0 it will move to "ct2". Rinse probe with DI water and then rinse with your second buffer.
6. Place in the second buffer. If you are doing a 2 pt., use the pH that is in your range (4 or 10).
7. Wait 30 seconds. Press [Enter] and wait for meter to show "ct3".
8. If doing a 2 pt. Calibration, press [M] twice to read samples.

Probe storage

Always store the probe in the solution of 3mol KCL solution. If not available, use pH 4.0

Contact us for supplies, equipment, and yearly certification options:

office@certify911.com

www.certify911.com